





# pCDH-CMV-MCS-EF1-copGFP-T2A-Puro

## 基本信息

出品公司	SBI
载体名称	pCDH-CMV-MCS-EF1-copGFP-T2A-Puro
质粒类型	慢病毒表达载体；cDNA 表达载体；双启动子载体
克隆方法	多克隆位点，限制性内切酶
启动子	CMV
载体大小	--
5' 测序引物及序列	CMV-F :CGCAAATGGGCGGTAGGCGTG
3' 测序引物及序列	--
载体标签	无
载体抗性	氨苄青霉素 (Ampicillin)
筛选标记	GFP、Puromycin
克隆菌株	E.coli cells(RecA-)推荐: Stbl3 ,OmniMAX 2 T1R
宿主细胞 (系)	常用细胞系, 如 HeLa, HEK293, HT1080, H1299
备注	pCDH-CMV-MCS-EF1-copGFP-T2A-Puro 慢病毒表达载体是基于 HIV 的慢病毒载体； 用于 cDNA 表达和克隆；高效转染细胞，建立稳定细胞系； CMV 启动子驱动目的基因的高水平表达，EF1a 启动子驱动报告基因的中等水平的表达。
产品目录号	CL525-01
稳定性	稳表达
组成型/诱导型	组成型
病毒/非病毒	慢病毒 (HIV)

## 载体质粒图谱和多克隆位点信息

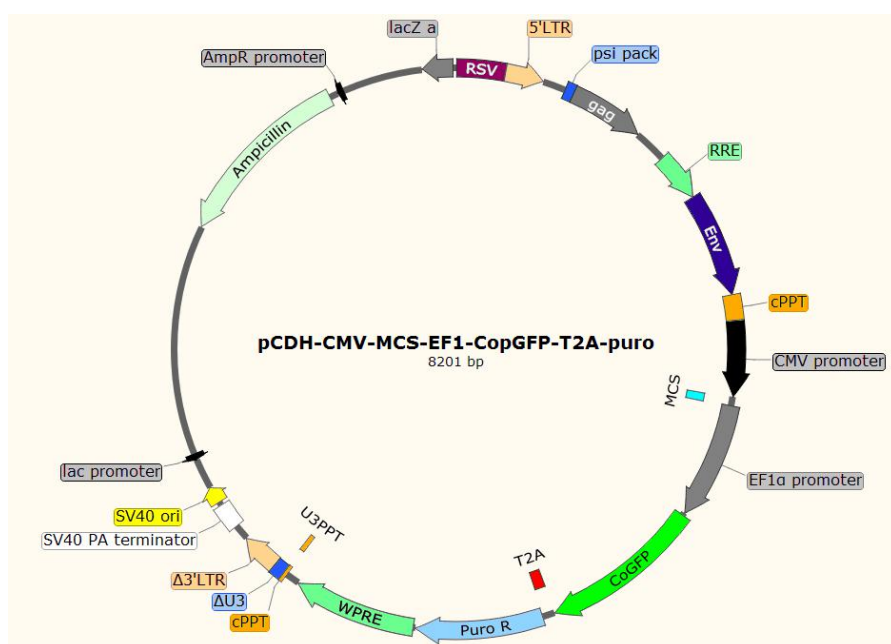

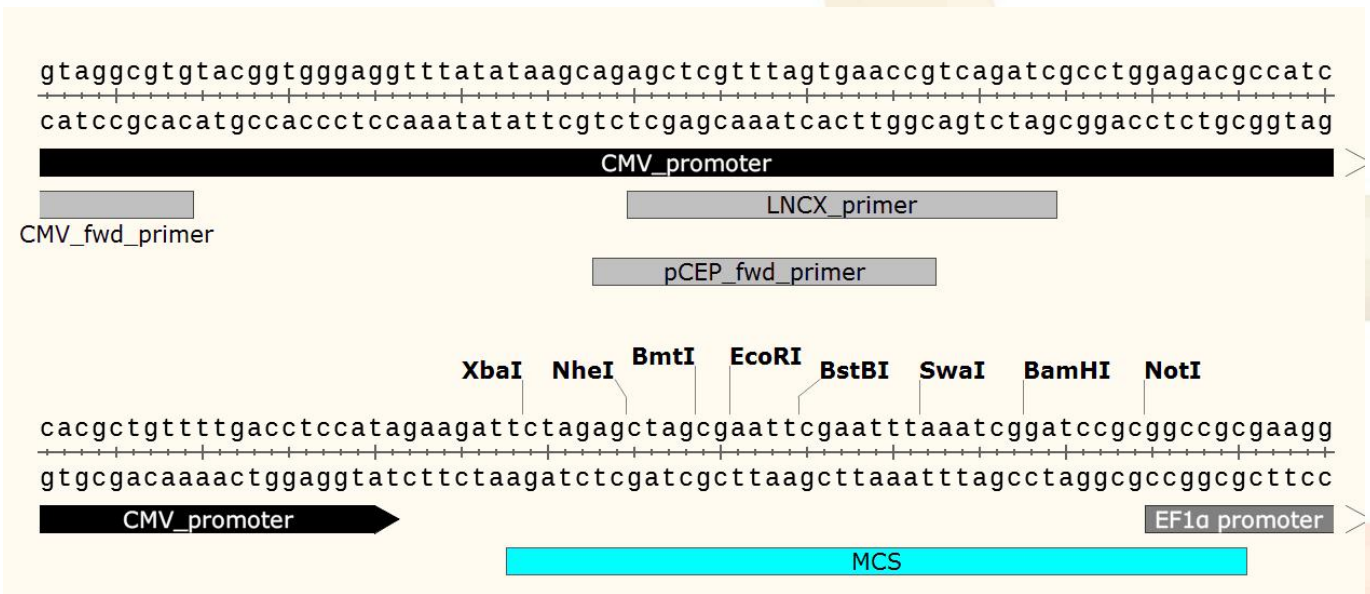

### 背景简介:

This manual provides details and information necessary to generate expression constructs of your gene of interest in the pCDH cDNA Cloning and Expression Lentivectors. Specifically, it provides critical instructions on amplification and cloning cDNA into the pCDH vectors, and verification of the final expression constructs. This manual does not include information on packaging the pCDH expression constructs into pseudotyped viral particles or transducing your target cells of choice with these particles. This information is available in the user manual Lentivector Expression Systems: Guide to Packaging and Transduction of Target Cells which is available on the SBI website. Before using the reagents and material supplied with this system, please read the entire manual.

### 基于 HIV-1 的 pCDH 慢病毒载体特征:

- Multiple Cloning Site (MCS)—for cloning the gene of interest in the MCS located downstream of the CMV promoter.
- WPRE element—enhances stability and translation of the CMV-driven transcripts.
- SV40 polyadenylation signal—enables efficient termination of transcription and processing of recombinant transcripts.
- Hybrid RSV/5LTR promoter—provides a high level of expression of the full-length viral transcript in producer 293 cells.
- Genetic elements (cPPT, gag, env, LTRs)—necessary for packaging, transducing, and stably integrating the viral expression construct into genomic DNA.
- SV40 origin—for stable propagation of the pCDH plasmid in mammalian cells.
- pUC origin—for high copy replication and maintenance of the plasmid in E.coli cells.
- Ampicillin resistance gene—for selection in E.coli cells.

### pCDH 慢病毒表达载体的优势:

Lentiviral expression vectors are the most effective vehicles for the delivery and expression of a gene of interest to almost any mammalian cell—including non-dividing cells and model organisms (C.A. Machida, 2003; M. Federico, 2003; W. C. Heiser, 2004). As with standard plasmid vectors, it is possible to introduce lentivector expression constructs in plasmid form into the cells with low-to-medium efficiency using conventional transfection protocols. However, by packaging the lentivector construct into viral particles, you can obtain highly efficient transduction of expression

constructs—even with the most difficult to transfect cells, such as primary, stem, and differentiated cells. The expression construct transduced in target cells is integrated into genomic DNA and provides stable, long-term expression of the target gene.

### **pCDH 慢病毒载体的包装载体及细胞系**

The expression lentivector contains the genetic elements responsible for packaging, transduction, stable integration of the viral expression construct into genomic DNA, and expression of the target gene sequence. The packaging vector provides all the proteins essential for transcription and packaging of an RNA copy of the expression construct into recombinant viral particles. To produce a high titer of viral particles, expression and packaging vectors are transiently co-transfected into producer mammalian cells (e.g., HEK 293 cells). For a detailed description of SBI's Lentivector expression system, please refer to the Lentivector Expression System user manual.

#### **启动子的选择:**

SBI provides a collection of cDNA cloning and expression vectors for various applications. A gene of interest can be cloned under a CMV or EF1 promoter with or without another expression cassette for a reporter gene (copGFP or PuroR). Genes can be either expressed transiently through transfection or stably expressed in a target cell line through transduction with packaged viral particles.

The major concern of cDNA expression in lentivectors is the efficiency level and stability of expression in target cell lines.

The Cytomegalovirus (CMV) promoter is a strong and most commonly used viral promoter that constitutively expresses downstream genes. While the CMV promoter works perfectly in the most common cell lines, it shows poor expression in some stem cell lines and hematopoietic cell lines (R.F. Doll, 1996; E.D. Papadakis, 2004).

The housekeeping elongation factor 1 $\alpha$  (EF1) promoter has been shown to exceed and outlast CMV-mediated expression in retroviral, lentiviral, and adenoviral vectors, in hematopoietic cell lines (K. Tokushige 1997; H. Nakai, 1998; C. Teschendorf, 2002). EF1 also performs well in most common cell lines.

MSCV promoter is the 5'-LTR promoter of murine stem cell virus. When a portion of the U3 region of the 3' HIV LTR was replaced with the U3 region of MSCV LTR, the resulted hybrid HIV/MSCV LTR has dramatically increased the transgene expression level in human CD34+ hematopoietic cells (J.K. Choi, 2001). After integration into genomic DNA, this promoter transcribes a long transcript with an intron in the 5'UTR flanked with splice donor and acceptor sites derived from the lentiviral vector. Further studies found that additional CpG mutations in the MSCV LTR reduced transcriptional silencing in embryonic stem cells (C.S. Swindle, 2004). We constructed cDNA expression vectors with the CpG-deficient MSCV incorporated into the 3' HIV LTR. After integration into genomic DNA, 3'MSCV/LTR will replace the 5'LTR and provide a high level of expression of the target gene and reporter gene downstream.

### **SBI 第三代慢病毒载体**

SBI offers a third generation of the most popular HIV-1 based lentivector expression system which consists of three main components:

- (1) The lentiviral expression vector (e.g., pCDH-EF1-MCS-T2A-Puro)
- (2) The lentiviral packaging plasmids (e.g., pPACKH1 Packaging Plasmid mix)
- (3) A pseudoviral particle producer cell line (e.g., 293TN cells)

#### **2A Peptide-enabled dual expression system**

Coexpression of a reporter gene together with a gene of interest is a useful approach for selecting transfected or transduced cells. This is commonly achieved by using two independent internal promoters, such as CMV and EF1 in pCDH-CMV-MCSEF1-copGFP, or by linking two transgenes with an internal ribosomal entry site (IRES) element in a single bicistronic transcript. Many dual promoter pairs have shown a high level of expression of both transgenes in standard cell lines— however, promoter interference often occurs in some cell lines. There are also two main problems that limit the use of IRES: the large size and the imbalanced expression between the first and second cistrons (H. Mizuguchi, 2000; X.Yu, 2003).

The “self-cleaving” 2A peptides have been used successfully to generate multiple proteins from a single promoter in many applications (P. de Felipe, 2004; M.J. Osborn, 2005; P. de Felipe, 2006). The 2A-like sequences exist in several viruses and are used to mediate protein cleavage from a single open reading frame. Through a ribosomal skip mechanism, the 2A peptide prevents normal peptide bond formation between the 2A glycine and the 2B proline without affecting the translation of 2B (M.L. Donnelly, 2001):SBI’s cDNA expression vectors incorporate the 2A-like sequence (T2A) from the insect virus *Thosea asigna* to mediate the coexpression of a reporter gene with the target cDNA. Reporter genes have been cloned at either the first or second positions, and we achieved high expression levels at both locations.

## 载体序列

```

1  acgcgtgtag tcttatgcaa tactcttgta gtcttgcaac atggtaacga tgagttagca
61  acatgcctta caaggagaga aaaagcaccg tgcatgccga ttggtggaag taagtggtga
121 cgatcgtgcc ttattaggaa ggcaacagac gggctctgaca tggattggac gaaccactga
181 attgccgcat tgcagagata ttgtatttaa gtgcctagct cgatacaata aacgggtctc
241 tctggttaga ccagatctga gcctgggagc tctctggeta actagggaac cactgctta
301 agcctcaata aagcttgcct tgagtgcttc aagtagtgtg tgcccgtctg ttgtgtgact
361 ctggtaacta gagatccctc agaccctttt agtcagtgtg gaaaatctct agcagtggcg
421 cccgaacagg gacctgaaag cgaaagggaa accagagctc tctcgacgca ggactcggct
481 tgctgaagcg cgcacggcaa gaggcgaggg gcggcgactg gtgagtacgc caaaaatfff
541 gactagcggg ggctagaagg agagagatgg gtgcgagagc gtcagtatta agcggggggg
601 aattagatcg cgatgggaaa aaattcgggt aaggccaggg ggaaagaaaa aatataaatt
661 aaaacatata gtatgggcaa gcagggagct agaacgattc gcagttaatc ctggcctggt
721 agaaacatca gaaggctgta gacaaatact gggacagcta caaccatccc ttcagacagg
781 atcagaagaa cttagatcat tatataatac agtagcaacc ctctattgtg tgcatcaaag
841 gatagagata aaagacacca aggaagcttt agacaagata gaggaagagc aaaacaaaag
901 taagaccacc gcacagcaag cggccactga tcttcagacc tggaggagga gatatgaggg
961 acaattggag aagtgaatta tataaatata aagtagtaaa aattgaacca ttaggagtag
1021 caccaccaa ggcaaagaga agagtgggtc agagagaaaa aagagcagtg ggaataggag
1081 ctttgttctt tgggttcttg ggagcagcag gaagcactat gggcgcagcc tcaatgacgc
1141 tgacggtaca ggccagacaa ttattgtctg gtatagtgca gcagcagaac aatttctga
1201 gggctattga ggcgcaacag catctgttgc aactcacagt ctggggcatc aagcagctcc
1261 aggcaagaat cctggctgtg gaaagatacc taaaggatca acagctcctg gggatttggg
1321 gttgctctgg aaaactcatt tgcaccactg ctgtgccttg gaatgctagt tggagtaata
1381 aatctctgga acagattgga atcacacgac ctggatggag tgggacagag aaattaacaa
1441 ttacacaagc ttaatacact ccttaattga agaatcgcaa aaccagcaag aaaagaatga
1501 acaagaatta ttggaattag ataaatgggc aagtttgtgg aattggttta acataacaaa
1561 ttggctgtgg tatataaaat tattcataat gatagtagga ggcttggtag gtttaagaat
1621 agtttttctg gtacttttcta tagtgaatag agttaggcag ggatattcac cattatcgtt
1681 tcagaccacac ctcccaaccc cgagggggacc cgacaggccc gaaggaatag aagaagaagg
1741 tggagagaga gacagagaca gatccattcg attagtgaac ggatctcgac ggtatcggtt
    
```



1801 aacttttaaa agaaaagggg ggattggggg gtacagtgca ggggaaagaa tagtagacat  
 1861 aatagcaaca gacatacaaa ctaaagaatt acaaaaacaa attacaaaat tcaaaatfff  
 1921 atcgatacta gtattatgcc cagtacatga ccttatggga ctttctact tggcagtaca  
 1981 tctacgtatt agtcatcgct attaccatgg tgatgcggtt ttggcagtac atcaatgggc  
 2041 gtggatagcg gtttgactca cggggatttc caagtctcca cccattgac gtcaatggga  
 2101 gtttgttttg gcaccaaaat caacgggact ttccaaaatg tcgtaacaac tccgccccat  
 2161 tgacgcaaat gggcggtagg cgtgtacggg gggaggttta tataagcaga gctcgtttag  
 2221 tgaaccgtca gatcgcctgg agacgccatc cacgctgttt tgacctccat agaagattct  
 2281 agagctagcg aattcgaatt taaatcggat ccgcggccgc gaaggatctg cgatcgtcc  
 2341 ggtgcccgtc agtgggcaga gcgcacatcg cccacagtcc ccgagaagt ggggggaggg  
 2401 gtcggcaatt gaacgggtgc ctagagaagg tggcgcgggg taaactggga aagtgatgtc  
 2461 gtgtactggc tccgcctttt tcccgagggt gggggagaac cgtatataag tgcagtatgc  
 2521 gccgtgaacg ttctttttcg caacgggttt gccgccagaa cacagctgaa gcttcgaggg  
 2581 gctcgcattc ctctttcacg cggccgccgc cctacctgag gccgccatcc acgccggtg  
 2641 agtcgcgttc tgccgcctcc cgctgtggt gcctctgaa ctgctccgc cgtctaggta  
 2701 agtttaaaagc tcaggctgag accgggcctt tgccggcgc tcccttgag cctacctaga  
 2761 ctacgccggc tctccacgct ttgcctgacc ctgcttgct aactctactg ctttgtttcg  
 2821 ttttctgttc tgcgccgta cagatccaag ctgtgaccgg cgctactgct agatggagag  
 2881 cgacgagagc ggctgcccc ccatggagat cgagtgccgc atcaccggca cctgaacgg  
 2941 cgtggagttc gagctgggtg gcggcggaga gggcaccccc aagcagggcc gcacgacaa  
 3001 caagatgaag agcaccaaag gcgccctgac ctccagcccc tactgtctga gccacgtgat  
 3061 gggctacggc ttctaccact tcggcaccta cccagcggc tacgagaacc ctttctgca  
 3121 cgccatcaac aacggcggct acaccaaac ccgcatcgag aagtacgagg acggcggcgt  
 3181 gctgcacgtg agcttcagct accgctacga ggccggccgc gtgatcggcg acttcaaggt  
 3241 ggtgggcacc ggcttcccc aggacagcgt gatcttcacc gacaagatca tccgagcaa  
 3301 cgccaccgtg gacacactgc accccatggg cgataacgtg ctggtgggca gcttcgccc  
 3361 caccttcagc ctgcgcgacg gcggctacta cagcttcgtg gtggacagcc acatgcactt  
 3421 caagagcgcc atccaccca gcacctgca gaacgggggc cccatgttcg ctttcgccc  
 3481 cgtggaggag ctgcacagca acaccgagct gggcatcgtg gattaccagc acgcttcaa  
 3541 gacccccatc gccttcgcca gatcccgcgc tcagtcgtcc aattctgccc tggacggcac  
 3601 cgccggacc ggctccaccg gatctcgtg agggcagagg aagtcttcta acatgcggtg  
 3661 acgtggagga gaatcccggc ctttccggga tgaccgagta caagcccag gtgcgcctcg  
 3721 ccaccgcga cgacgtcccc agggccgtac gcacctcgc cgccgcttc gccactacc  
 3781 ccgccacgcg ccacaccgtc gatccggacc gccacatcga gcgggtcacc gagctgcaag  
 3841 aactcttctt cacgcgcgtc gggctcgaca tcggcaaggt gtgggtcgcg gacgacggcg  
 3901 ccgcggtggc ggtctggacc acgccggaga gcgtcgaagc gggggcgggtg ttcgccgaga  
 3961 tcggcccgcg catggccgag ttgagcgggt cccggctggc cgcgagcaa cagatggaag  
 4021 gcctctggc gccgcaccgg cccaaggagc ccgcgtggtt cctggcccacc gtcggcgtct  
 4081 cgcccacca ccagggaag ggtctgggca gcgccgtct gctccccgga gtggagcggg  
 4141 ccgagcgcgc cgggggtgcc gccttcttgg agacctccgc gccccgaac ctccccctt  
 4201 acgagcggct cggttcacc gtcaccgcg acgtcaggt gccgaagga ccgcgcacct  
 4261 ggtgcatgac ccgcaagccc ggtgcctgag tcgacaatca acctctggat taaaaattt  
 4321 gtgaaagatt gactggtatt cttaactatg ttgctcttt tacgtatgt ggatacgtg  
 4381 ctttaatgcc tttgtatcat gctattgctt cccgtatggc tttcattttc tctcttgg  
 4441 ataaatcctg gttgctgtct ctttatgagg agttgtggcc cgttgtcagg caacgtggcg  
 4501 tggtgtgcac tgtgtttgct gacgcaacc ccactggtg gggcattgcc accacctgtc  
 4561 agctcctttc cgggactttc gctttcccc tccctattgc cacggcggaa ctatcggc  
 4621 cctgccttgc ccgctgctgg acaggggctc ggctgttggg cactgacaat tccgtggtg

4681 tgtcggggaa atcatcgtcc tttccttggc tgctcgctg tgttgccacc tggattctgc  
4741 gcgggacgtc cttctgctac gtcccttcgg cctcaatcc agcggacctt ccttcccgcg  
4801 gcctgctgcc ggctctgcgg cctcttccgc gtcttcgcct tcgccctcag acgagtcgga  
4861 tctccctttg ggccgcctcc ccgcttgta cctttaagac caatgactta caagcgagct  
4921 gtagatctta gccacttttt aaaagaaaag gggggactgg aagggtaat tcaactccaa  
4981 cgaaaataag atctgctttt tgcttgtaact gggctctctt ggtagacca gatctgagcc  
5041 tgggagctct ctggctaact agggaaacca ctgcttaagc ctcaataaag cttgccttga  
5101 gtgcttcaag tagtgtgtgc ccgtctgttg tigtactctg gtaactagag atccctcaga  
5161 cccttttagt cagtgtggaa aatctctagc agtagtagtt catgtcatct tattattcag  
5221 tatttataac ttgcaaagaa atgaatatca gagagtgaga ggaacttggt tattgcagct  
5281 tataatgggt acaataaag caatagcatc acaaatttca caaataaagc atttttttca  
5341 ctgcattcta gttgtggttt gtccaaactc atcaatgat cttatcatgt ctggctctag  
5401 ctatcccgcc cctaactccg cccagttccg cccattctcc gccccatggc tgactaattt  
5461 tttttattta tgcagaggcc gaggccgcct cggcctctga gctattccag aagtagtgag  
5521 gaggtttttt tggaggccta gacttttgca gagacggccc aaattcgtaa tcatggctat  
5581 agctgtttcc tgtgtgaaat tgttatccgc tcacaattcc acacaacata cgagccggaa  
5641 gcataaagtg taaagcctgg ggtgcctaata gtagtgagta actcacatta attgcgttgc  
5701 gctcactgcc cgctttccag tcgggaaacc tgcctgcca gctgcattaa tgaatcgccc  
5761 aacgcgcggg gagaggcggg ttgcgtattg ggcgctcttc cgcttctcgc cteactgact  
5821 cgctgcgctc ggtcgttcgg ctgcggcgag cggatcagc tcaactcaaag gcgtaatac  
5881 ggttatccac agaatcaggg gataacgcag gaaagaacat gtgagcaaaa gcccagcaaa  
5941 aggccaggaa ccgtaaaaag gccgcgttgc tggcgttttt ccataggctc cgccccctg  
6001 acgagcatca caaaaatcga cgctcaagtc agaggtggcg aaaccgaca ggactataaa  
6061 gataccaggc gtttccccct ggaagctccc tcgtgcgctc tctgtttccg accctgccc  
6121 ttaccggata cctgtccgcc tttctccctt cgggaagcgt ggcgctttct catagctcac  
6181 gctgtaggta tctcagttcg gtgtaggtcg ttcgctccaa gctgggctgt gtgcacgaac  
6241 cccccgttca gcccgaccgc tgcgccttat ccgtaacta tcgtcttgag tccaaccg  
6301 taagacacga cttatcgcca ctggcagcag ccactggtaa caggattagc agagcgaggt  
6361 atgtaggcgg tgctacagag ttcttgaagt ggtggcctaa ctacggctac actagaagga  
6421 cagtatttgg tatctgcgct ctgctgaagc cagttacctt cggaaaaaga gttgtagct  
6481 cttgatccgg caaacaacc accgctggta gcggtggttt ttttgttgc aagcagcaga  
6541 ttacgcgcag aaaaaaagga tctcaagaag atcctttgat cttttctacg ggtctgacg  
6601 ctcagtggaa cgaaaactca cgtaaggga ttttggatcat gagattatca aaaaggatct  
6661 tcacctagat ctttttaaat taaaaatgaa gttttaaact aatctaaagt atatatgagt  
6721 aaacttggtc tgacagttac caatgcttaa tcagttaggc acctatctca gcgatctgtc  
6781 tatttcgttc atccatagtt gcctgactcc ccgtcgtgta gataactacg atacgggagg  
6841 gcttaccatc tggccccagt gctgcaatga taccgcgaga cccacgctca ccggtccag  
6901 atttatcagc aataaaccag ccagccggaa gggccgagcg cagaagtggc cctgcaactt  
6961 tatccgcctc catccagtct attaattggt gccgggaagc tagagtaagt agttgccag  
7021 ttaatagttt gcgcaacggt gttgccattg ctacaggcat cgtggtgtca cgctcgtcgt  
7081 ttggtatggc ttcatcagc tccggttccc aacgatcaag gcgagttaca tgatccccc  
7141 tgttgtgcaa aaaagcgggt agctccttcg gtctccgat cgttgtcaga agtaagtgg  
7201 ccgcagtgtt atcactcatg gttatggcag cactgcataa ttctcttact gteatgceat  
7261 ccgtaagatg cttttctgtg actggtgagt actcaaccaa gtcattctga gaatagtga  
7321 tgcggcgacc gagttgctct tgccccgct caatacggga taataccgcg ccacatagca  
7381 gaactttaaa agtgctcatc attggaaaac gttcttcggg gcgaaaactc tcaaggatct  
7441 taccgctggt gagatccagt tcgatgtaac ccactcgtgc acccaactga tcttcagcat  
7501 cttttacttt caccagcgtt tctgggtgag caaaaacagg aaggcaaat gcccaaaaa

7561 agggaataag ggcgacacgg aatgttgaa tactcatact cttcettttt caatattatt  
7621 gaagcattta tcagggttat tgtctcatga gcggatacat atttgaatgt atttagaaaa  
7681 ataaacaaat aggggttccg cgcacatttc cccgaaaagt gccacctgac gtctaagaaa  
7741 ccattattat catgacatta acctataaaa ataggcgtat cacgaggccc tttcgtctcg  
7801 cgcgtttcgg tgatgacggt gaaaacctct gacacatgca gctcccggag acggtcacag  
7861 cttgtctgta agcggatgcc gggagcagac aagcccgtca gggcgcgtca gcgggtgttg  
7921 gcgggtgtcg gggctggctt aactatgcgg catcagagca gattgtactg agagtgcacc  
7981 atatgcggtg tgaaataccg cacagatgcg taaggagaaa ataccgcac aggcgccatt  
8041 cgccattcag gctgcgcaac tgttgggaag ggcgatcggg gcgggcctct tcgctattac  
8101 gccagctggc gaaaggggga tgtgctgcaa ggcgattaag ttgggtaacg ccagggtttt  
8161 ccagtcacg acgttgtaaa acgacggcca gtgccaagct g

BM20211208